

# Computer-Controllable High-Current Universal LED Controllers

(Part Numbers: SLC-HA02-U, SLC-HV02-U)

## PRODUCT DESCRIPTION

### FEATURES

- ◆ Computer controllable
- ◆ Universal - suitable for any LED
- ◆ Capable of driving variable loads
- ◆ User friendly application software with GUI
- ◆ SDK for custom applications
- ◆ Normal, Strobe and Trigger mode for every channel
- ◆ Programmable constant current, pulse-width modulation and/or arbitrary waveform
- ◆ Driving current up to **2000mA** in DC mode and up to **3500mA** in pulse mode, with over current protection
- ◆ Up to 11.5V output voltage for each channel
- ◆ Programmable rising or falling edge external trigger
- ◆ Built-in non-volatile memory, can be used without a PC

### APPLICATIONS

- ◆ Machine vision
- ◆ Displays
- ◆ Microscopy
- ◆ Semiconductor equipment
- ◆ Testing instruments
- ◆ Medical instruments
- ◆ Lighting

Mightex has developed a series of computer-controllable, multi-channel, universal LED controllers, which can be used to drive any type of LED in any of the three (3) modes: 'NORMAL' (or 'constant current'), 'STROBE', and/or external 'TRIGGER' mode. Each unit comes with PC-based software with a user-friendly GUI, which enables users to drive LEDs without the need to write any code. In addition, a powerful SDK is provided, in order for users to write their own software and to integrate Mightex's LED drivers into their own systems. Furthermore, the drivers have a built-in security feature, allowing users to limit LED driving current and voltage.



This datasheet covers product series SLC-HA and -HV, which currently include 2 models (i.e. SLC-HA02-U, SLC-HV02-U). The following table, which can also be used as a product selection guide, compares the key features of the 2 product models.

P/N	# of Channels	Control Mode <sup>(1)</sup>			Arbitrary Waveform <sup>(2)</sup>	Interface	Forward Voltage Monitoring
		NORMAL	STROBE	TRIGGER			
SLC-HA02-U	2	•	•	•		USB	
SLC-HV02-U	2	•	•	•		USB	•

Notes: (1) Each output channel can be individually configured to work in one of the following three (3) modes, controlled through a PC-based software with GUI. In all three modes, overdrive current limit can be set:

**Normal:** Constant current output at any value from 0mA to 2000mA with 1mA resolution.

**Trigger:** External trigger signal could be used to turn on each individual channel, generating driving current with any user-defined waveform. Alternatively, each output channel can work under the "FOLLOWER" mode, in which the current output follows the waveform of the trigger input; and

**Strobe:** Internal Strobe Generator generates frequencies as high as 25KHz. The strobe signal (i.e. current levels, duty cycle and strobe frequency) can be set through software.

# Computer-Controllable High-Current Universal LED Controllers

(Part Numbers: SLC-HA02-U, SLC-HV02-U)

## PERFORMANCE SPECIFICATIONS

### ELECTRICAL SPECIFICATION:

Parameters	SLC-HAxx-x	SLC-HVxx-x	Unit
Power Supply Input Voltage, V(dc)	9 ~ 12		V
Power Supply Input Current	4,000		mA
Per Channel Driving Voltage (max)	Up to 11.5		V
Maximum Per Channel Driving Current	2,000 ("NORMAL" MODE)		mA
	3,500 ("STROBE" or "TRIGGER" MODE)		mA
Output Current Resolution	1		mA
Output Current Linearity	+/-4 (or +/-0.5%)		mA
Output Current Repeatability	+/-1 (or +/-0.2%)		mA
Trigger Input High Level	3.3 ~ 10.0		V
Trigger Input Low Level	0.8 (Max.)		V
Forward Voltage Monitoring Accuracy	N/A	+/-10	mV

### TIMING SPECIFICATION:

Parameters	SLC-HAxx-x	SLC-HVxx-x	Unit
Timing Resolution	20		µs
# of Data Points for Waveform Definition	2		
Trigger Pulse Width	100 (Minimum)		µs
Max Trigger Delay	25		µs


## OPERATION CONDITION

Operating Temperature Range: 0°C ~ 45°C  
 Storage Temperature Range: -25°C ~ 85°C  
 Relative Humidity, Non-condensing: 5% ~ 95%

## DIMENSION AND WEIGHT

Dimension: 201mm(L) x 147mm (W) x 40mm (H)  
 Weight: 600g


## INSTALLATION DRAWING

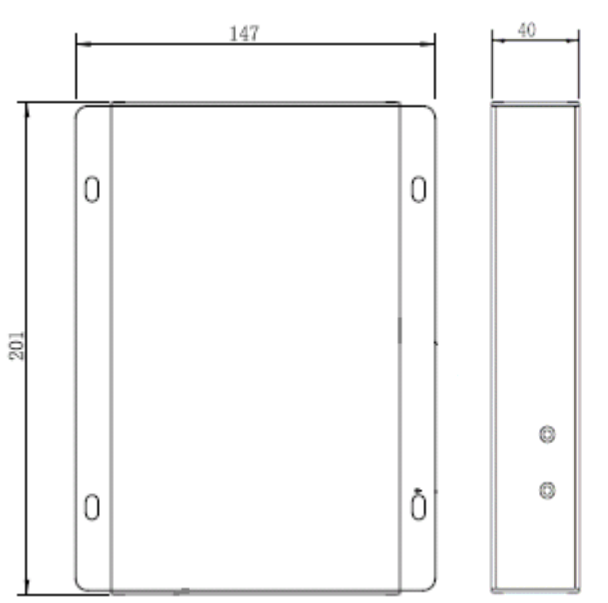

# Computer-Controllable High-Current Universal LED Controllers


(Part Numbers: SLC-HA02-U, SLC-HV02-U)

## APPLICATION DIAGRAM

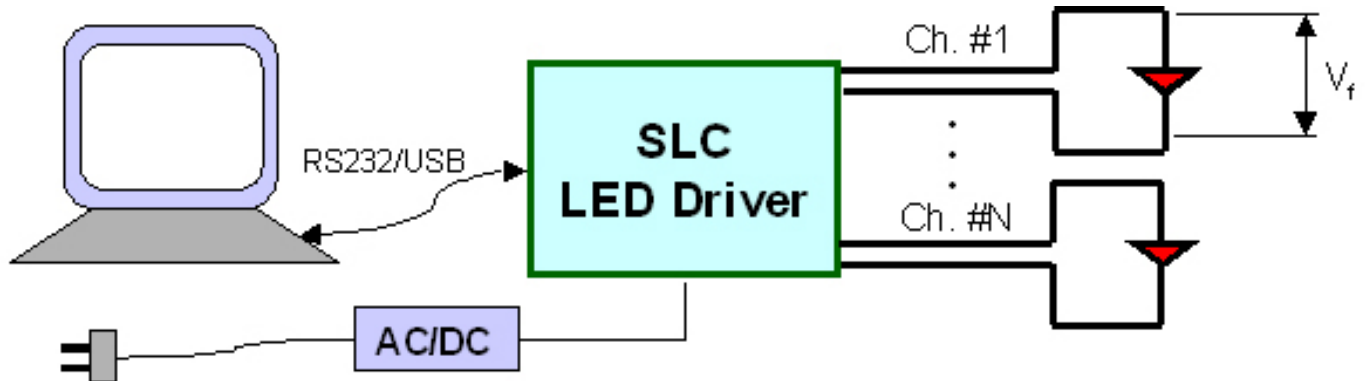

## EXAMPLE OF GRAPHICAL USER INTERFACE

- GUI application for manual channel control
- Pre-programmed patterns for Trigger and Strobe mode

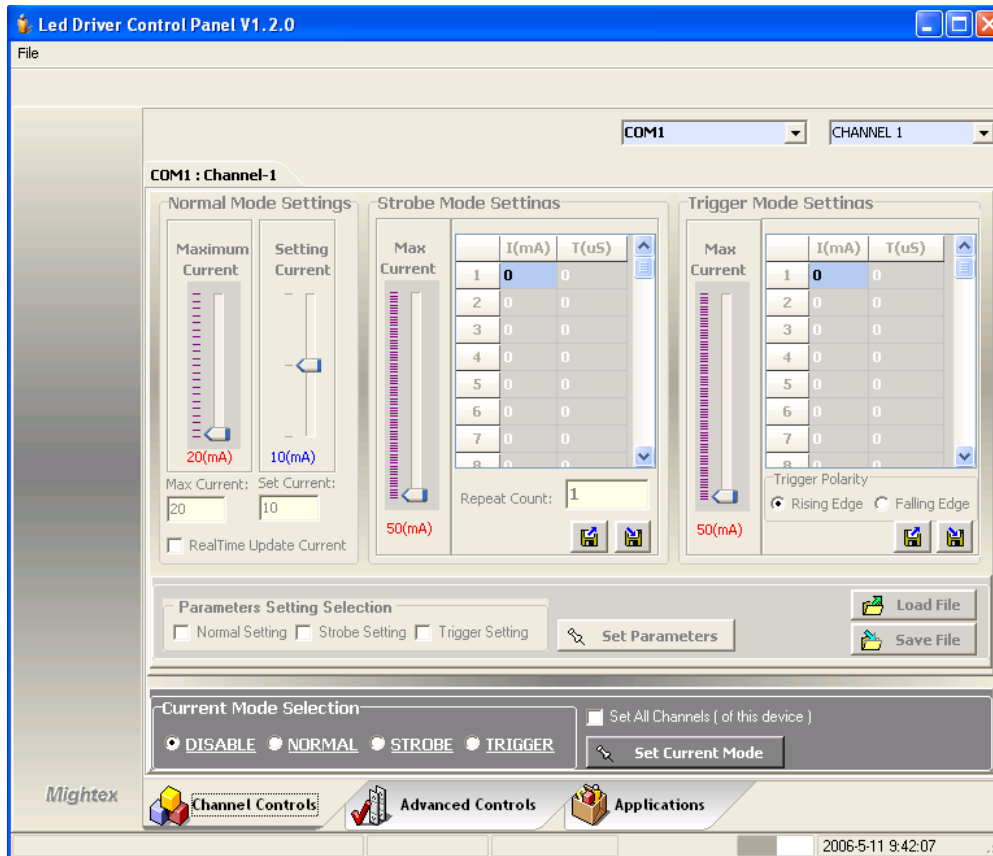

For customized solutions, please call 1-416-840 4991 or email [sales@mightex.com](mailto:sales@mightex.com).